

## Updating the firmware 340 prepASH

### Downloading and unpacking

1. Download the actual firmware package, available in the download section of our website <http://www.precisa.com>.  
The package comes as zip folder named "340prepASH\_Firmware-Package\_Vx-x-x.zip", e.g. "340prepASH\_Firmware-Package\_V1-0-4.zip".
2. For unpacking just doubleclick the downloaded zip folder and let the operating system, Windows 7 or higher, do this automatically. **Important: Do NOT unzip by means of "WinZip" or similar applications; the files could be damaged and unusable!**

The unzipped folder consists of:

prepASH\_updater: Folder with file "signature.txt" and Zip-File "prepASH\_updater.zip".

Installation instructions: This file, named "Installation\_Instruction.pdf".

Release notes: Text file named "ReleaseNotes\_Vx-x-x.pdf" listing all release notes about this firmware version.

### Updating

1. Copy the folder "prepASH\_updater" onto an empty USB stick.
2. Insert the USB stick into the prepASH terminal.
3. Login with a user that has the necessary permissions to update the prepASH.
4. Navigate to the menu <Settings><About device><Software> and press the <Update> button.
5. Once the process is complete, the application will restart automatically.