

## Release Notes 1.3.3

### New features

### Miscellaneous

- Remote command list included in context help of setting "Data exchange – ..."
- Soft button "1/...d" renamed to "x Decimals"

### Bug fixes

- Info item "ID": Text-initialisation when changing application

## Release Notes 1.3.2

### New features

- Application "Dynamic" add-on "Checkweigher" included
- Soft button "1/...d" for resolution-reduction included

### Miscellaneous

- Add-on "Peak Hold" protocol item "Peak value" included
- Add-on "Statistics" action button "Max. n" included
- MSW setting "Capture below minimum" included
- For Statistics- and Protocol-Load change adjustable "Min. Load" included
- Data exchange mode "Measurement transfer" transfer item "ID" included
- Remote control commands "D<text>", "DN" and "@<text>", "@N" for writing text into the display included  
(Remote command list is now included in context help "Data exchange – Serial Port")

### Bug fixes

- French text corrected

## Release Notes 1.3.1

### New features

- Printer LP4024, printout in Chinese language available (no settings on printer required).  
Note: Printer Firmware LP 4024 Vers:3.0 required

# NEWS



Max 225g Min 1mg d=0.01/0.1mg e=1mg

Administrator - 22.05.2024 16:02

## Miscellaneous

- Translations revised
- Context help for addon Identification added

## Bug fixes

- Add-On Identification: Customizing of soft buttons [ID1..6]
- Add-On Checkweigher: Function of soft button [Capture nominal]

## Release Notes 1.3.0

### New features

- Time synchronisation via the network added, works from terminal version BSP V2.2.4 upwards (see device information).
- Time zone setting added. Automatic daylight-saving time regulation included.

### Miscellaneous

- Data output (serial port, mass storage and network) in Chinese language encoded as character set GB18030 for traditional Chinese and GB2312 (subset of GB18030) for simplified Chinese.

## Bug fixes

## Release Notes 1.2.4

### New features

### Miscellaneous

## Bug fixes

- RTC deviation correction

## Release Notes 1.2.3

### New features

### Miscellaneous

- Service remote command CLRHISTORYcccc added.
- Public remote command FWHISPRINT added.

### Bug fixes

## Release Notes 1.2.1

### New features

### Miscellaneous

- Writing to a CSV file revised. CSV separator selectable, semicolon or comma

### Bug fixes